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The screenshot displays the software's procedure assignment window. On the left, a list of 'Procedures Available' includes 1ML.SPE, 3ML example.SPE, 3ML.SPE, CLEAN.SPE, Flow Test slow speed.spe, Flow Test.spe, Load 0,5mL SLE+.spe, Load 0,5mL SLE+.spe, Load 1,1mL SLE+.spe, Load 1mL SLE+.spe, and Purge and clean.spe. On the right, 'Procedures assigned to:' shows 'Module 0' with samples 1 through 10. Below this is a 'Delete Selected Items' button. At the bottom, there are checkboxes for Module 0 through 9, all of which are checked, along with 'Check All' and 'Uncheck All' buttons. A 'Close' button is at the bottom left, and 'OK / Save' and 'Help' buttons are at the bottom right.

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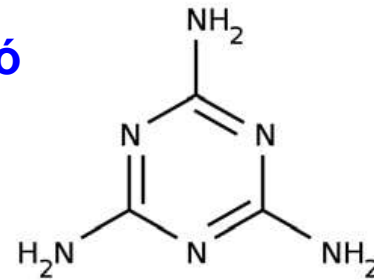
* Biotage Application Note AN749

Extraction of Melamine from Milk (Infant Formula) using EVOLUTE® CX Columns and RapidTrace+ Workstation

BIOTAGE RAPIDTRACE+

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

	Step	Source	Destination	Volume (mL)	Flow (mL/min)
1	Condition	Methanol	Organic Waste	3.0	2.00
2	Condition	Ammonium acetate	Organic Waste	3.0	2.00
3	Load	Sample	Biological Waste	2.0	1.00
4	Purge-Cannula	Ammonium acetate	Organic Waste	2.0	30.00
5	Rinse	Ammonium acetate	Organic Waste	3.0	2.00
6	Rinse	Methanol	Organic Waste	3.0	2.00
7	Collect	Methanol:NH ₄ OH (95:5)	Fraction 1	3.0	1.00
8	Purge-Cannula	Ammonium acetate	Organic Waste	6.0	30.00
9	Purge-Cannula	Methanol	Cannula Waste	6.0	30.00
10	Purge-Cannula	Water	Cannula Waste	2.0	30.00

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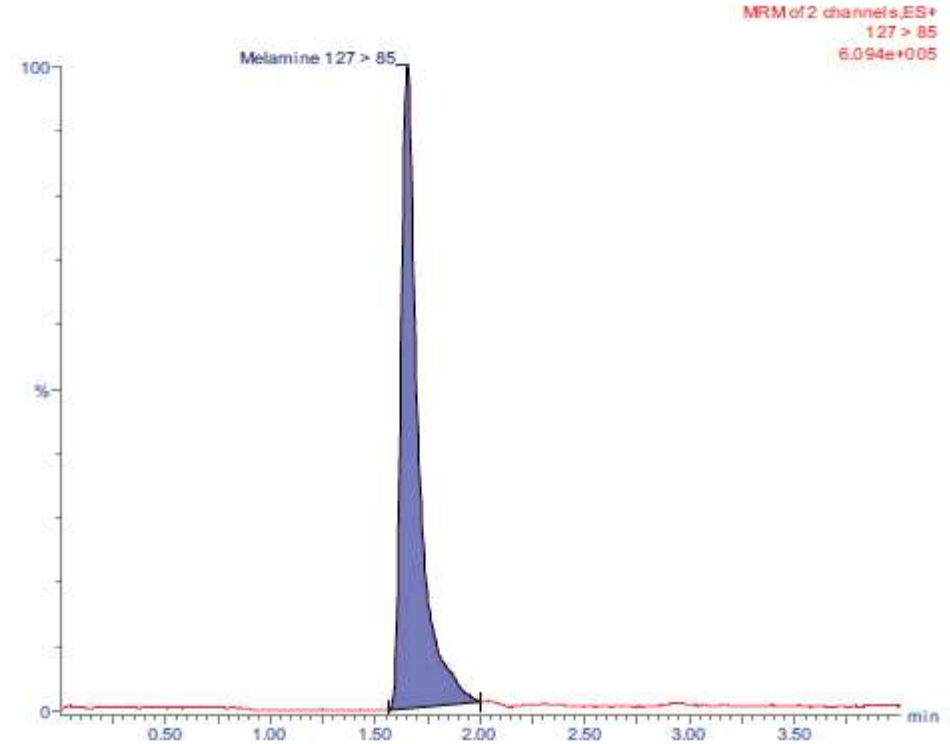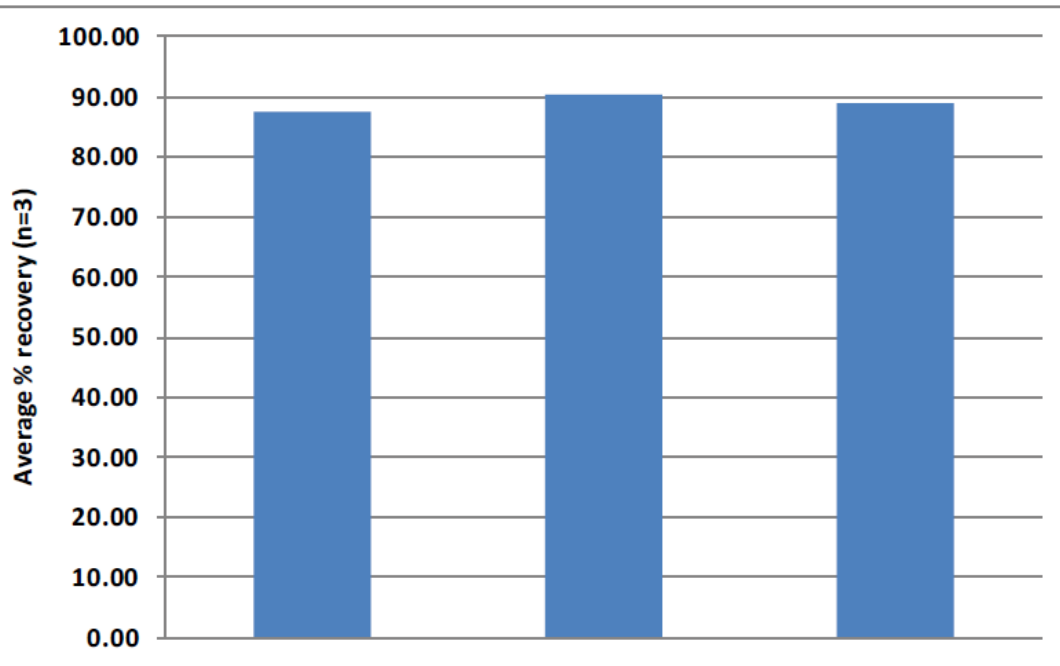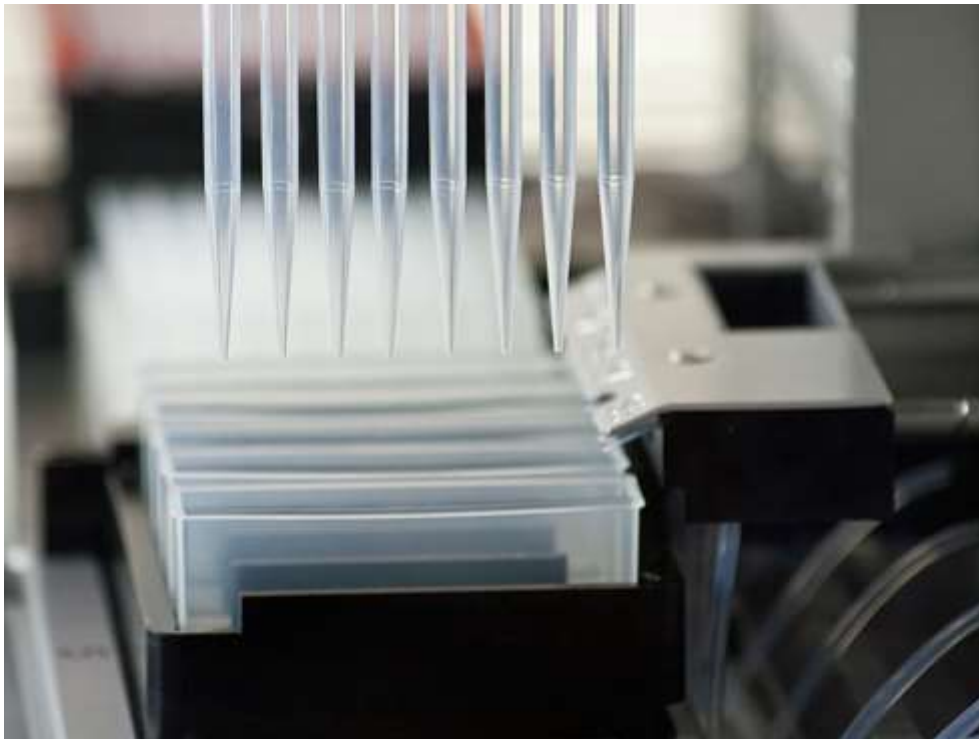
BIOTAGE EXTRAHERA

< Back View SPE Method - EVOLUTE ABN (non polar), ...

Method name: EVOLUTE ABN (non polar), 6 mL columns

Sample plate/rack: 12 x 75 mm test tubes, 24

Extraction media: 6 mL column rack, 24

Pretreatment: On

Conditioning: Off

Equilibration: Off

Sample Load: On

Wash: On

Elution: On

Number of steps: 1 Pressure (bar): 1,5 Plate dry after last elution? Yes Plate dry time (s): 120 Dispose tips after each step? No

Solvent: Methanol

Volume (µL): 2500

Positive pressure time (s): 120 Advanced..

Repeat (number of times): 2 Pause after this step? No

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< Back Prepare Run - EVOLUTE CX EXPRESS Run >

Sample Plate/Rack ID: Extraction Media Lot No.: Collection Plate/Rack ID:

Solvent Tips: 1

Extraction Media: 3

Solvent Reservoirs: 5

Sample Tips: 2

Sample Plate/Rack: 4

Carousels: 6

Report and Comment:

Carousal Setup:

Extraction Media:

Extraction Media:	Extraction plate, 96
Sample Plate/Rack:	1 mL sample plate, 96
Solvent Tip Columns Needed:	3
Sample Tip Columns Needed:	0
Total Solvent Consumption:	
0.03M AcOH, pH 4	0.000 mL
Methanol	0.000 mL
MeOH-NH4OH (95:5)	0.000 mL
Pretreatment:	
0.03M AcOH, pH 4	750 µL
Conditioning:	
	Not used
Equilibration:	
	Not used
Sample Load:	
Aqueous sample	500 µL
Wash:	
0.03M AcOH, pH 4	1000 µL
Methanol	1000 µL
Elution:	
MeOH-NH4OH (95:5)	1000 µL

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* Biotage Product Note PPS358, Product Note PPS360
Biotage® Extrahera™ Precision and accuracy Performance

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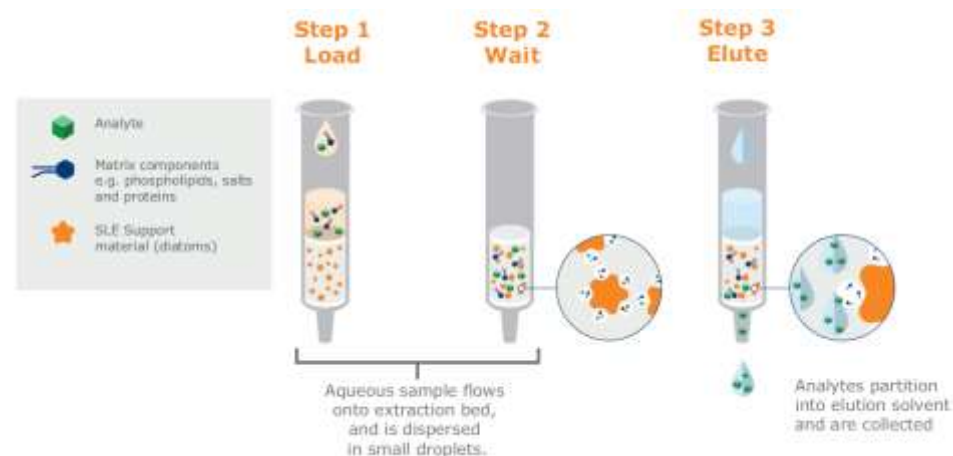
* Biotage poster presented at TIAFT, Brisbane, 2016
Evaluation of Novel Automated Sample Preparation Compared to Manual Processing in
Forensic Toxicology

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